



Formaldehyde Spill Response Kit Directions

Kit Number: (48625, 48630)

Kit Contents:

- Pair of Nitrile Gloves
- Respirator
- Open Back Gown
- Vented Goggles
- FSC-1 Solidifier (11oz. Bottle)
- Scoop/Scraper (Quantity: 4)
- Wiper Pads (Quantity: 3)
- Clear Disposal Bags (Quantity: 4 / Dimensions: 30" x 36")
- Twist Ties (Quantity: 4)
- and Instructions

Directions:

- 1.) Put on open back gown, respirator, vented goggles and gloves.
*Always wear protective gear and exercise caution during clean up.
- 2.) Sprinkle FSC-1 over spill evenly. The fluid will quickly set; add more FSC-1 as necessary.
- 3.) Remove solidified material with scoop and scraper and place in one of the supplied white disposal bags
- 4.) Clean area with a hard surface wipe or disinfectant (**For best results use Safetec's® SaniZide Plus® surface disinfectant spray or wipe**). Wipe up any residue left behind with supplied wiper pads.
- 5.) After procedure is completed remove gloves, vented goggles, respirator and open back gown....Place all contaminated materials in white disposal bag. Seal bag with supplied twist tie.
- 6.) Dispose of all waste in accordance with local, state, and federal regulations.

SAFETEC OF AMERICA, Inc. Toll Free: 1-800-456-7077

Address: 887 Kensington Ave. Buffalo, NY 14215 Web: www.safetec.com